

# **Instructions for Plasma Cleaner:**

Before starting, always make sure the sample is loaded correctly. The compartment must be unhooked and placed over the sample. Be careful not to crack the glass on the metal base.

## **STEPS:**

- 1) Open door, lift the lid if not up, and place the sample inside.
- 2) Turn on the Vacuum (green power switch to the left of the plasma cleaner)
- 3) Turn on System Vacuum (above the cleaner)
- 4) Turn on Gas Control
- 5) Adjust dial for SCFH Air until it is around 3
- 6) Run plasma cleaner for 3 minutes
- 7) Turn off Gas Control, System Vacuum, and Vacuum
- 8) Unload Sample and leave Glass Compartment on the hook.